


Rotational movement:  
 ~30 arc-sec/  
 10 degrees of adjustment of YAW.

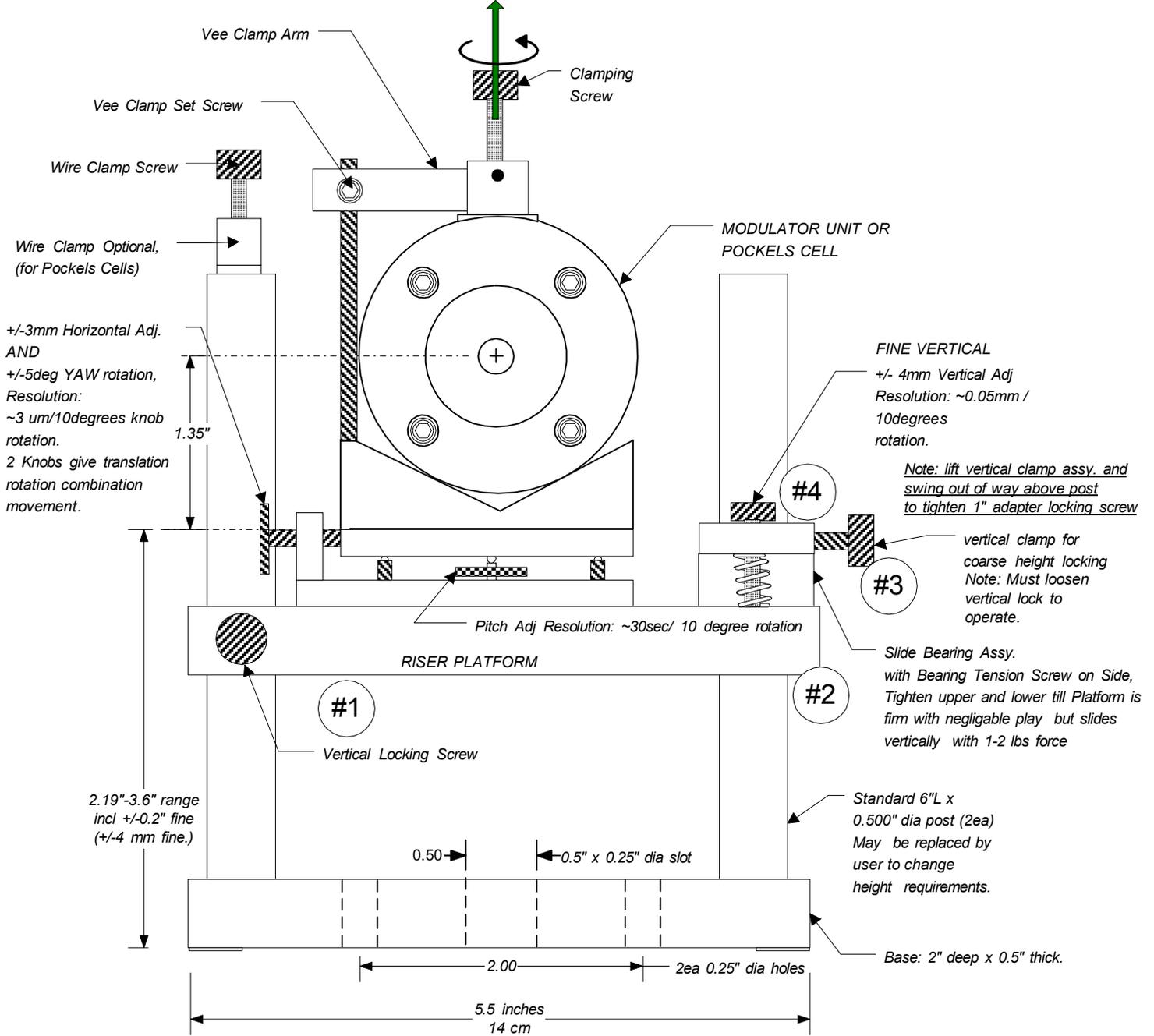

- A. To adjust Coarse Vertical, loosen #1.
- B. Loosen #3
- C. Lift or Lower Riser Platform by gripping, #2, till aperture is at or close to beam height
- D. Lock #3
- E. Adjust fine height, #4, till set
- F. Lock height in place, #1

- A. To adjust Fine Vertical, loosen #1.
- B. Adjust fine height, #4, till set
- C. Lock height in place, #1

QUANTUMT TECHNOLOGY, INC. LAKE MARY, FLA, USA 32746			
MODEL: MTM MOUNT WITH MODULATOR			
SIZE	FSCM NO	DWG NO	REV
A		MODEL-MTM-MOUNT WITH MODULATOR	C
SCALE	1 : 1	SHEET	1 OF 1